



The ERWEKA

Export Manager

for direct PDF and XML export



ERWEKA

Overview

Introducing the ERWEKA Export Manager: Simple and Fast Data Export! With the ERWEKA Export Manager, users can seamlessly export test protocols from devices SVM II, TAR II, EasyCheck, MultiCheck 6 and ZT 720/30 via their network, effortlessly converting them into PDF and XML formats. The Windows software, once installed on any PC within your lab network, automatically transforms your test report data into PDF and XML files. Say goodbye to time-consuming manual scanning of printed reports.

But that's not all. Our user-friendly interface allows you to define your desired destination folder, whether it is a local drive or a designated network location. Integrated into the labs own IT network, security measures and data compliance protocols valid for the network are also valid for those folders, ensuring valuable data always remains safeguarded. No compromises on data security or integrity.

Each exported file is specifically named, combining the device type, serial number, and the date and time of data reception. This means users will never lose track of valuable data, maintaining an effortless overview for easy accessibility and analysis.

The ERWEKA Export Manager is a light-weight solution, catering to both new and existing equipment. With its remarkable ability to receive data from multiple devices simultaneously, users will save time, effort, and money and streamline their laboratory workflows.

Supported devices

The Export Manager software supports the following ERWEKA devices:

- ZT 72x (from Firmware-Version 1.20)
- ZT 73x (from Firmware-Version 2.02)
- EasyCheck (from Firmware-Version 1.38)
- MultiCheck 6 (from Firmware-Version 1.38)
- SVM II (from Firmware-Version 1.15)
- TAR II (from Firmware-Version 1.01)

Status

Shows the status of the network connection.

- Red: Values for "Port number" and "Target folder" are missing or incorrect.
- Yellow: The values entered are correct.
- Green: Data reception has been started. The software is ready to receive data from the device.

IP address

Shows the IP address of the PC on which the Export Manager is running. It is determined automatically.

Port number

Sets the number of the port where data is received from the device.

Supported data formats

The software supports ASCII and XML data formats for receiving data, and PDF and XML data formats for exporting.



Quick and easy
PDF / XML export



Numerous
supported
ERWEKA devices



USB stick with
software & copy
protection dongle

Start listening

Switches the Export Manager to the "Waiting" status, in which data can be received. The button is active only after you have entered correct values for "Port number" and "Target folder".

File Help Export Manager

ERWEKA

Status Stop listening

IP address 185.210.***

Port number 85**

Target folder C:\Users\Erweka\Desktop ...

Event screen

[Server] waiting for a client...

[Server] Client connected: /172.22.0.100:12850

ERWEKA Device: ZT 722 Page 1/1

Direct measurement report

OQ status: invalid PV date: 01.05.2022
Device S/N: 134728 Firmware version: 01.19.0

Test started: 07.02.2022 17:07
Test ended: 07.02.2022 17:22

Batch: Machine:
Stability: Info:
Basket number: Basket type: Type A

Test station: 1
Test status: Passed
Pre-Set temperature: 37.5°C
Nominal speed: 30 DPM
Pre-Set runtime: 00:30:00
Total runtime: 00:13:56

Min-Max Temperature: 37.1-37.6°C Min-Max Speed: 29.7-30.5 DPM

Vial	Result
1	00:11:46
2	00:09:37
3	00:10:11
4	00:13:53
5	00:13:56
6	00:08:40

Statistics
Number of passed samples: 6
Minimal time: 00:08:40 Maximal time: 00:13:56
Average time: 00:11:20 Max - Min: 00:05:16
Std.Deviation: 00:02:13 Rel.Deviation: 19.6%

Test result: Passed

Performed by: ERWEKA Reviewed by: _____
ERWEKA Staff Date: _____

Copyright © Erweka Version 1.00.0

Target folder

Sets the folder path where the protocol will be saved. Click to set a path.

Event screen

Displays information about the status of the Export Manager and the received data (live, for control only).



Contact

Are you curious and want to find out more?
Head over to our website and download our product brochures,
watch videos of our equipment in action or find the ERWEKA
dealer of your country.



E-Mail: sales@erweka.com
Phone: +49 6103 92426-200
Fax: +49 6103 92426-999



support@erweka.com



www.erweka.com



spareparts@erweka.com



www.facebook.com/erweka.gmbh

ERWEKA GmbH

Pittlerstr. 45
63225 Langen
Germany

E-Mail: sales@erweka.com
Phone: +49 6103 92426-200
Fax: +49 6103 92426-999

Technical specifications of products described are
stated without warranty and subject to change at
any time without further notice. **v.1.0.6.23**

ERWEKA